

# ULT25NEU Quick Guide:



Download Operator and Quick Guide manuals in the language of choice at: [www.stirlingultracold.com/manuals](http://www.stirlingultracold.com/manuals)


-----clip, laminate, keep with freezer-----

## INITIAL SETUP

Port Stopper/Sealant	<b>!</b>	The Port Stopper/Sealant must be in place during operation.
Power Indicator	<b>A</b>	The unit is ON when lit
Power ON	<b>B</b>	Press and hold for two beeps
Power OFF	<b>B</b>	Press and hold for three beeps
LCD Display	<b>C</b>	Displays chamber temperature
Display Setpoint	<b>F</b>	Press/Release to display current setpoint, S ## displayed
Return to Temp Display	<b>F</b>	Press/Release or wait eight seconds
Change Setpoint	<b>F</b>	Press and hold until <i>Set T LED</i> (E) glows, S ## displayed, then ▲ or ▼ (D)

## ADVANCED FUNCTIONS

Change to Preset	Press and hold <i>Set</i> (F) until <i>Set T LED</i> (E) glows, S ## displayed, then select <i>Preset 1, 2 or 3</i> (J)
Postpone Over Temp Alarm	When <i>Over Temp LED</i> (I) glows, press <i>Preset 3</i> (J) until time shows (1.0 h) (If <i>Over Temp LED</i> is not glowing, then <i>Preset 3</i> acts as a Temp Preset)
Return to Temp Display	Wait eight seconds, press <i>Preset 3</i> to return to temperature display
See Error Code Again	If <i>Error LED</i> (H) is glowing, press <i>Preset 2</i> (J) briefly (If <i>Error LED</i> is not glowing, then <i>Preset 2</i> becomes Temp Preset)
Lock Panel	Press and hold <i>Preset 1 AND 2 AND 3</i> (J) until <i>Locked LED</i> (G) turns on (~five seconds)
Unlock Panel	Press and hold <i>Preset 1 AND 2 AND 3</i> (J) until <i>Locked LED</i> (G) turns off (~five seconds)
LCD Display Codes	P ## – Preset Temp, ## h – Alarm Silenced, E ## – Error Code (see operating manual)

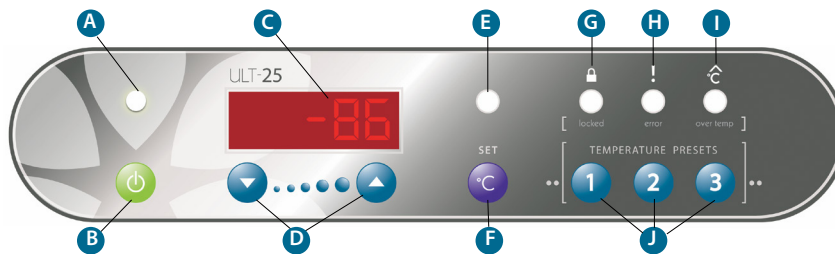

- When unit is first turned on, red Over T light will be on until temperature is within 10° of set point.
- Panel cannot be locked until temperature is within 10° of set point.



6000 Poston Road Athens, Ohio 45701, USA t: 740.274.7900 / 855.274.7900 f: 740.274.7901 [BioLifeSolutions.com](http://BioLifeSolutions.com)

©2023 Stirling Ultracold, a part of BioLife Solutions. All Rights Reserved.  
Global Cooling technology is manufactured under U.S. and International patents.  
Stirling Ultracold is a trademark of Global Cooling, Inc.  
Specifications subject to change without notice. Refer to [BioLifeSolutions.com](http://BioLifeSolutions.com) for the latest specifications.

